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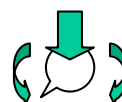
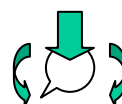
Please use after reserves.

Call 7472 and tell the use time. (Sorry,Japanese only)

How to use of centrifuge model 7780 II.

1. Adjust the balance of weight of sample. Set it in the rotor symmetrically.
2. Turn on the power switch(The left side of main body) of the centrifuge.
3. Open the door with the OPEN button.
4. Set the rotor on the axis. The cap of rotor is closed.
(Do again if the cap of rotor doesn't close.)
5. The door is closed.
6. Press NORMAL button.
7. Turn the dial, and set the parameter. (TEMP,SPEED and TIME)
The “temp limit” lamp lights when TEMP dial is pressed.
Centrifuge stops at preset temperature +5℃ when the “temp limit” lamp lights.
If SPEED dial is pressed, it is converted from “rpm” to “g”. (or opposite)
The unit of time changes if TIME button is pressed.
8. Set acceleration and deceleration with ACCEL,DECEL button. The rapid acceleration/deceleration when a lot of numbers of LED lights.
9. Press the START button to start centrifuge.
10. It stops automatically when time comes. Take out the rotor.
(Turn off power with the door opened.)
11. Write the record in the notebook.

Press a dial

Turn a dial

Wash the rotor in distilled water when the sample spills in the rotor.